


A Simple Setup to use Semi-Preparative HPLC Columns on Biotage® Selekt

The newly released Biotage® Selekt flash chromatography instrument can be run at a maximum flowrate of 300 mL/min or a maximum pressure of 30 bar. These high flowrates and pressures enable a user to perform chromatography using not only dry-packed, single-use plastic flash columns containing small ($\geq 20 \mu\text{m}$) spherical silica particles, but also semi-preparative, slurry-packed HPLC columns for multiple use with smaller ($\leq 20 \mu\text{m}$) spherical silica particles.

To use semi preparative columns on Selekt you can build your own kit, and a simple way can be by using the list below (most of the components available at www.IDEX-hs.com).

Injection Valve, 2 Position-6-Port .063 – Red

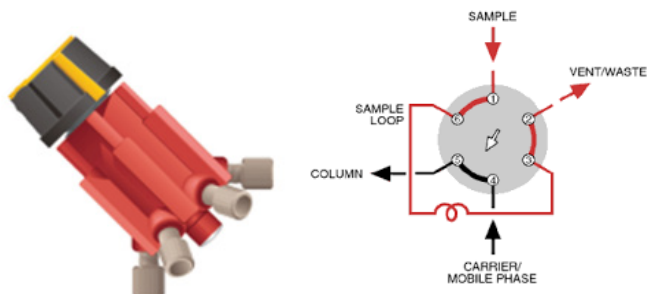

Super Flangeless Fittings



**Loop 3 mL Steel 1/8" OD, 0.080" ID
with Super Flangeless Nuts**



Sample Inject Port
Luer Adapter Female Luer to 1/4-28 Male, Delrin®



**Column Tubing 1/16" OD Steel, 0.030" ID with
HPLC Connectors and Super Flangeless Connectors**



EUROPE

Main Office: +46 18 565900
Toll Free: +800 18 565710
Fax: +46 18 591922
Order Tel: +46 18 565710
Order Fax: +46 18 565705
order@biotage.com
Support Tel: +46 18 56 59 11
Support Fax: +46 18 56 57 11
eu-1-pointsupport@biotage.com

NORTH & LATIN AMERICA

Main Office: +1 704 654 4900
Toll Free: +1 800 446 4752
Fax: +1 704 654 4917
Order Tel: +1 704 654 4900
Order Fax: +1 434 296 8217
ordermailbox@biotage.com
Support Tel: +1 800 446 4752
Outside US: +1 704 654 4900
us-1-pointsupport@biotage.com

JAPAN

Tel: +81 3 5627 3123
Fax: +81 3 5627 3121
jp_order@biotage.com
jp-1-pointsupport@biotage.com

CHINA

Tel: +86 21 68162810
Fax: +86 21 68162829
cn_order@biotage.com
cn-1-pointsupport@biotage.com

KOREA

Tel: +82 31 706 8500
Fax: +82 31 706 8510
korea_info@biotage.com
kr-1-pointsupport@biotage.com

INDIA

Tel: +91 22 4005 3712
india@biotage.com

Distributors in other regions
are listed on www.biotage.com

Literature Number: UI425

© 2019 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. A list of all trademarks owned by Biotage AB is available at www.biotage.com/legal. Other product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.