

Optical Rotation Measurement for High-concentration Acidic Sample Solutions

< Key Points >

- The P-4000 polarimeters offer high durability for acidic sample measurements because of an acid-resistant coating on the sample compartment wall.

< Introduction >

In certain optical rotation tests described in the United States pharmacopeia (USP) and the Japanese pharmacopoeia (JP), sample preparation requires the use of acidic solutions such as 5 N hydrochloric acid (USP) or 6 mol/L hydrochloric acid (JP). Because optical rotation is temperature-dependent, it is important to monitor the temperature during measurements. The P-4000 series polarimeters are equipped with a temperature sensor that can be inserted directly into the cell. This cell sensor has an acid-resistant coating, allowing direct measurement of the temperature of acidic sample solutions. To protect the instrument when measuring acidic liquid samples, the P-4000 series incorporates the following features:

1. Acid-resistant coating on the sample compartment wall
2. Quartz window preventing acidic gas from entering the instrument*
3. Ventilation of acidic gas*
4. Connection to a ventilation duct**

* Standard feature only for the P-4200 and P-4300

** Optional for the P-4200 and P-4300

This application note describes the measurement of L-alanine in 6 mol/L hydrochloric acid.

< Keywords >

Polarimeter, acid resistant, 6 mol/L hydrochloric acid, L-alanine



Fig. 1 P-4200 Polarimeter

< Experimental >

● Sample

L-alanine (research reagent)

In accordance with the monograph described in JP XVIII¹⁾ (after drying, 2.5 g, 6 mol/L hydrochloric acid TS, 25 mL, 100 mm), the sample was prepared by dissolving dried L-alanine in a 6 mol/L hydrochloric acid solution.

● System

Instrument: P-4200 polarimeter

Cell: RQC-450 8.5-mm rectangular quartz cell (path length: 100 mm)

● Measurement conditions

Light source: Na lamp Wavelength: D-line

Integration time: 5 sec Number of repeats: 1

< Results and Discussion >

Table 1 shows the measurement results for the specific rotation of L-alanine. The acid-resistant cell sensor enables accurate monitoring of the sample temperature, even for a strong acid solution of 6 mol/L hydrochloric acid, allowing precise measurement of the optical rotation.

Table 1 Results for L-alanine

Specific rotation	Temperature during measurement / °C
+14.824	20.4

< Conclusions >

The P-4000 series polarimeters allow accurate measurement of optical rotation for highly acidic sample solutions, such as 6 mol/L hydrochloric acid, while directly monitoring the sample temperature. Additionally, these polarimeters exhibit high resistance to acidic samples, enabling reliable operation while maintaining the instrument in optimal condition.

< References >

1. Ministry of Health, Labour and Welfare: June 7, 2021, MHLW Ministerial Notification No. 220, “The Japanese Pharmacopoeia 18th edition”, (2021).

Applicative Solution Lab Division Y. Kondo

This document has been prepared based on information available at the time of publication and is subject to revision without notice. Although the contents are checked with the utmost care, we do not guarantee their accuracy or completeness. JASCO Corporation assumes no responsibility or liability for any loss or damage incurred as a result of the use of any information contained in this document. The company names, product names, service marks and logos mentioned herein are the trademarks and registered trademarks of the respective companies. Copyright and other intellectual property rights in this document remain the property of JASCO Corporation. Please do not attempt to copy, modify, redistribute, or sell etc. in whole or in part without prior written permission.